

WARNING NOTICE: The experiments described in these materials are potentially hazardous and require a high level of safety training, special facilities and equipment, and supervision by appropriate individuals. You bear the sole responsibility, liability, and risk for the implementation of such safety procedures and measures. MIT shall have no responsibility, liability, or risk for the content or implementation of any of the material presented. [Legal Notices](#)

## ELECTROCOMPETENT *RHODOCOCCUS ERYTHROPOLIS* AN12

---

### Competent cells for electroporation

1. (Day or two before) Inoculate small (5-10 ml) LB or NBYE cultures of *Rhodococcus erythropolis* AN12; grow at 30°C
2. Transfer 0.1-5 ml of overnight culture of *Rhodococcus* to 200 ml NBYE, 0.05% Tween-80 in a 1L baffled flask
3. Incubate shaking at 30°C overnight or until O.D.<sub>600</sub> is between 0.8 and 1.0
4. Pellet the cells by centrifuging for 5 min at 6 000 rpm in a GSA rotor using sterile centrifuge bottles or 50 ml conical tubes and proper adapters (may have to spin twice to pool)
5. Resuspend the cell pellet in 30 ml ice-cold distilled, sterile water; transfer cells to sterile 50 ml conical tube (if you haven't already done so); Recentrifuge as in step 4 and discard the supernatant
6. Resuspend cell pellet one more time in distilled, sterile water; centrifuge as before; discard supernatant
7. Wash pellet once in 10 ml ice-cold, sterile 10% glycerol; centrifuge as before except at 8000 rpm; discard supernatant
8. Resuspend final cell pellet in 1 ml or less of ice-cold, sterile 10% glycerol
9. Aliquot 150 µl into sterile microfuge tubes and store at -80°C

### Electroporation of *Rhodococcus*

1. Thaw aliquots of electrocompetent *Rhodococcus* cells on ice
2. Mix DNA with 70µl cells in a sterile microfuge tube and incubate on ice for 5 min.
3. Electroporate DNA at 2.5 kV, 25 µF and 400 Ω
4. Immediately add 300 µl NBYE
5. Incubate cells for recovery at 30°C for 1-6 hours
6. Spread cells onto plates with appropriate antibiotics

#### NBYE

0.8% nutrient broth (Difco)

0.5% yeast extract (Difco)

#### Tween-80 stock

5ml Tween-80

95 ml distilled water

Filter sterilize